

SHORT GUIDE TO INSTALLING SPECTRAGENIUS

- Installation Instructions -

**Please follow the installation instructions as explained in this guide.
The whole process takes only a few minutes to complete.**

CONTENT

-	<i>INSTALLATION INSTRUCTIONS –</i>	1
1.	<i>GENERAL INFORMATION</i>	3
1.1	Time limited demo	3
1.2	Demo databases included	3
2.	<i>BEFORE INSTALLATION</i>	3
3.	<i>INSTALLATION</i>	3
3.1	Starting Installation	3
3.2	Installation of SpectraGenius	4
3.3	Starting SpectraGenius	6
4.	<i>UNINSTALLING SPECTRAGENIUS</i>	7
5.	<i>WHAT TO DO WHEN YOU HAVE QUESTIONS</i>	8

1. General Information

This installation guide describes how to install the SpectraGenius demo version from S.T. Japan-Europe GmbH. The S.T. Japan-Europe installer (Installer or Setup) endeavors to make the installation of S.T. Japan-Europe spectra databases as transparent and decision free as possible. Please follow the steps to install S.T. Japan-Europe spectra databases as described in the following sections.

1.1 Time limited demo

The SpectraGenius demo version will work for 30 days after installation. It is limited to saving 200 files.

1.2 Demo databases included

SpectraGenius comes with two demo databases in Thermo Galactic format. Each includes 200 ATR or Raman spectra, respectively.

2. Before Installation

- SpectraGenius will work only on 64-bit Windows 8 and higher versions.
- SpectraGenius will not work in a virtual machine (e.g., VMWare).
- Make sure that all your current work is saved properly and close all programs before beginning the installation.
- If you are using an Anti-Virus software, like e.g. Microsoft's Anti-Virus, Norton Anti-virus, or Central Point's Anti-Virus, it is recommended that you turn the program off before installing S.T.Japan-Europe SpectraGenius. You can turn it back on after installation.

3. Installation

3.1 Starting Installation

After preparing the system as described in the “Before Installation” section, unzip the SpectraGenius demo installer onto the PC/Laptop on which you want to install the spectra databases. Then, double click the **SpectraGenius_install_55.exe** file.

The *Welcome* dialog is displayed.

The *Cancel* button will stop the installer and will allow you to terminate the installation process without adding or changing any files. You can click the *Cancel* button at any time of the installation process.


The *Next* button continues with the installation of the spectra databases and the *Options* dialog is displayed.



The *Options* dialog shows a check box list with the following options:

- **Install SpectraGenius** – Installs SpectraGenius, the demo databases, and some demo spectra
- **Service** – Opens a Service Dialog

Please **do not** check the *Service* box. The *Service* box only needs to be checked if and when you are advised by the S.T.Japan-Europe Support Team to do so.

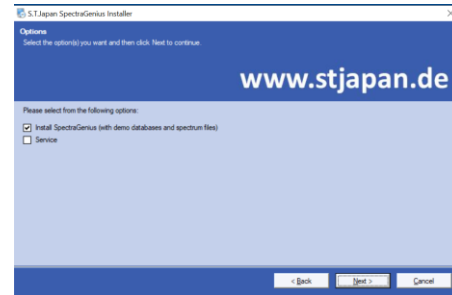

3.2 Installation of SpectraGenius

After clicking the *Next* button on the Options dialog, the installation of SpectraGenius starts, and the *License Agreement* is displayed.

Please note: If you wish to terminate the installation process at any time, click the *Cancel* button. This will stop the installer and allow you to terminate the installation process without adding or changing any further files.

When the *Back* button is selected, the installer will back up one installation step and display the previous installation dialog. The installation process continues by selecting the *Next* button.

If you disagree with terms and conditions of the License Agreement, click the *Cancel* button. This will stop the installer and allow you to terminate the installation process without adding or changing any files on your system.

If you agree with the Terms and Conditions of the License Agreement, click in the circle next to “*I agree to the terms of this license agreement*”, this will activate the *Next* button. To continue with the installation of SpectraGenius, click on the *Next* button. The installer displays the *Installation Folder* dialog. We recommend using the default install location. Click *Next*.

The progress bar indicates the status of the SpectraGenius installation. As the software is installed, the progress bar will indicate which file is currently being installed until all files are completely installed.

The installer for the demo databases will launch automatically. A couple of advertisement screens will appear – click *Next* on each one until you reach the license agreement for the demo databases.

If you disagree with terms and conditions of the License Agreement, click the *Cancel* button. This will stop the installer and allow you to terminate the installation process without adding or changing any files on your system.

If you agree with the Terms and Conditions of the License Agreement, click in the circle next to “*I agree to the terms of this license agreement*”, this will activate the *Next* button. To continue with the installation of the spectra databases click on the *Next* button.

On the next screen, please fill in your name and organization to receive more information on S.T. Japan-Europe’s spectra databases, then click *Next*.

The installer displays the *Installation Folder* dialog.

IMPORTANT: For optimum support, it is strongly recommended that the spectra databases are installed into the default installation folder C:\SpecLibs, which later will be referred to as *Database Installation Folder*. This folder will be created automatically during the spectra database installation. However, it is also possible to select another folder during the installation of the spectra databases.

To continue with the installation of the spectra databases, click the *Next* button.




The installer displays the *Ready to Install* dialog.

Continue the installation by selecting the *Next* button. The installer will display the *Installing* dialog.

The installation of the databases begins, and the databases are installed into the previously selected destination folder.

The progress bar indicates the status of the database installation. As the databases are installed, the progress bar will indicate which database is currently being installed until all purchased databases are completely installed.

If you wish to terminate the installation process during this process, click the *Cancel* button at any time. This will stop the installation process and allows you to terminate the installation process.

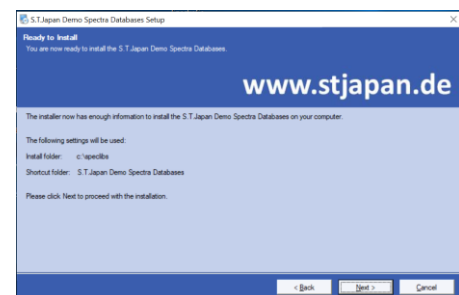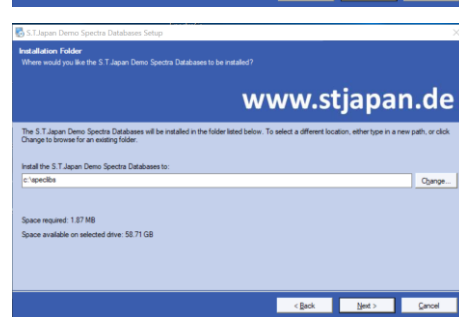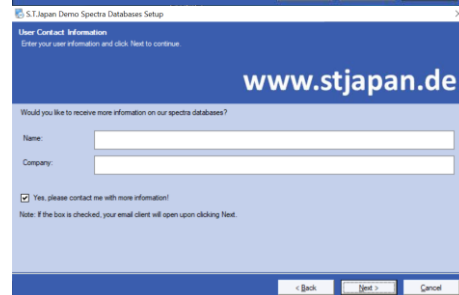


After the installation has been completed, the installer will display the *Installation Successful* dialog.

The spectra databases have been completely installed. To complete the installation process, click the *Finish* button. The installer will terminate, and the demo spectra installer will launch.



The installer displays the *Installation Folder* dialog. We recommend using the default path.


Click *Next* to continue.

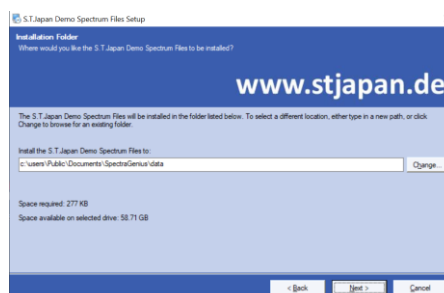

The installer displays the *Ready to Install* dialog.

Continue the installation by selecting the *Next* button. The installer will display the *Installing* dialog.

The installation of the demo files begins, and the files are installed into the previously selected destination folder.

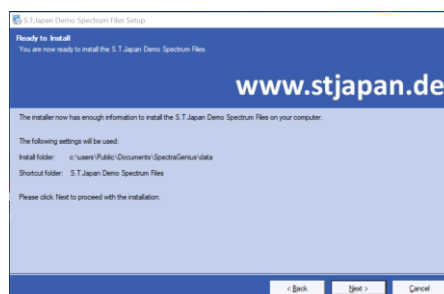


After the installation has been completed, the installer will display the *Installation Successful* dialog.

Everything has been completely installed. To complete the installation process, click the *Finish* button. The installer will terminate, and the Windows desktop will show.




3.3 Starting SpectraGenius


Clicking the SpectraGenius desktop icon will launch the SpectraGenius software.

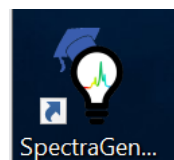


To access the Demo databases, click *Library Search* in the lower left area of the SpectraGenius screen, then click *Select Libraries*.

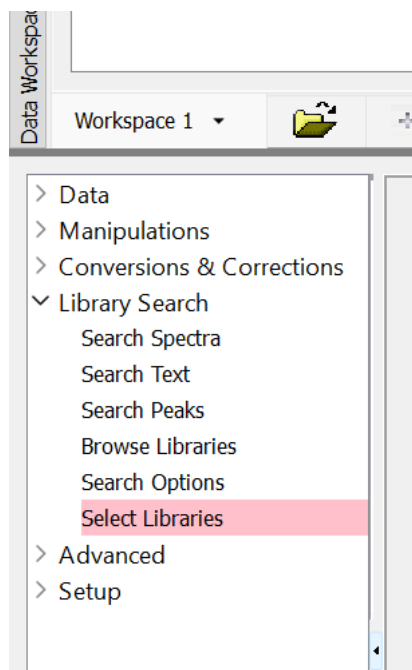


Under “All available libraries”, click *Add Library to Available List...* and then select *Add SpectralID library*. A file selection box will open. In it, browse to the folder where the demo databases have been installed (by default: c:\speclibs)

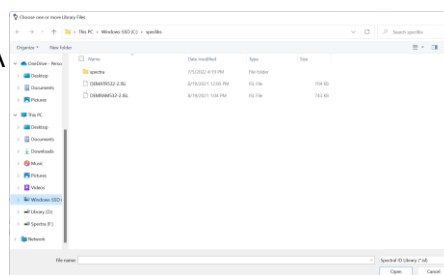

Select the one or both DEM*.ISL files, then click *Open*. The selected file(s) will appear in the list under “All available libraries”.

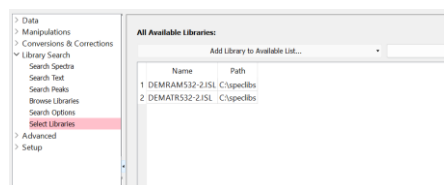

Select both entries, then click *Add Libraries to Search List* to make them available for use. The selected entries will appear in the list on the right.

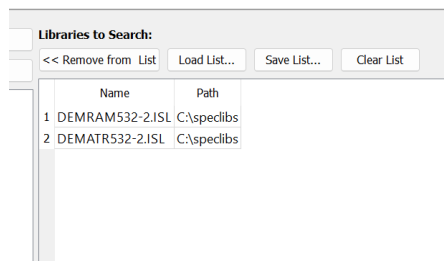

Done! For instructions on how to perform a search, or browse a library, please refer to the SpectraGenius help.

4. Uninstalling SpectraGenius

There are different ways to uninstall the spectra databases.

- **Uninstall from the Windows Start menu:** Select Start and look for the SpectraGenius Installer in the list shown. Then press and hold (or right-click) on SpectraGenius Installer and select Uninstall.

- **Uninstall from the Settings page:** Select Start, then select Settings > Apps > Apps & features. Select the SpectraGenius Installer, and then select Uninstall.
- **Uninstall from the Control Panel (for programs):** In the Windows search box on the taskbar, type Control Panel and select it from the results. Select Programs > Programs and Features. Press and hold (or right-click) on SpectraGenius Installer and select Uninstall or Uninstall/Change. Then follow the directions on the screen.

5. What to do when you have questions

If you have questions or experience difficulties with the installation of spectra databases, please do not hesitate to contact us for assistance

support@stjapan.de or call us at +49 (0)2234 956372.